

## DT8034

# 8 Channel 6 kV/ 1 mA Desktop Power Supply Module (USB/Ethernet/Touchscreen)



## Features

- 8 independent channels in a Desktop case (110/220V AC Powered)
- 6 kV / 1 mA output range
- Channels with either positive or negative polarity
- SHV coaxial output connectors
- Low Ripple
- Under/over-voltage alert, overcurrent and max. voltage protection
- Interlock logic for board enable and Individual channel kill
- 5 nA Current monitor resolution (with x10 Imon-Zoom: 500 pA)
- 2.8" color touch screen display
- Local and Remote control (USB2.0/Ethernet)
- Software Tools for easy channel management

## Description

The CAEN **Mod. DT8034** provides **8** independent 6 kV/ 1 mA channels in a desktop package (110/220V AC Powered). Each channel can provide a **6 kV** max voltage and a **1 mA** max current. The unit is available with positive, negative and "mixed" (4 positive and 4 negative) channels.

Channels outputs are delivered through SHV connectors. The HV output RAMP-UP and RAMP-DOWN rates may be selected independently for each channel in the range 1÷500 V/s in 1 V/s steps. The module features 5 nA Iset resolution.

### SHV connector

HV coaxial connector, 8 channels for Mod. DT8034

Consult our **connectors reference page** for technical information.

Module control can take place either **locally** thanks to a **2.8" Touchscreen Graphic color LCD display** with a completely redesigned user interface or **remotely**, via **USB 2.0** or **Ethernet**. A complete set of **Software Tools** is available to control these units; the user can freely download low level libraries, LabVIEW driver and Graphical application software.

Safety features allows the module to perform as a current generator and includes:

### Overvoltage and Undervoltage warning

When the output voltage differs from the programmed value by more than 2% of set value (minimum 10V).

### Overcurrent detection

When a channel attempts to exceed the programmed current limit, it signalled to be in "overcurrent" and enter in a TRIP status. The output voltage is varied to keep the current below the programmed limit for a programmable TRIP time, then the channel is switched off. If TRIP is set to "constant current mode", the channel behaves like a current generator.

### Hardware VMAX and IMAX

Maximum output voltage and maximum current value can be fixed, via front panel potentiometer, at the same common value for all the board channels. IMAX and VMAX values can be read out via software.

### Safety Board Interlock

Common Interlock logic for channels enable/disable and individual inputs signal for channel Kill function.

## Technical Specifications

### Package

Desktop module housed in a 239x84x184 mm<sup>3</sup> (WxHxD) Aluminium case  
(110/220V AC Powered)

### Output channels

8 channels, SHV connector  
Positive, Negative or Mixed (4-8 positive and 4-8 negative) Polarity; common ground

### Output Voltage

0 ÷ 6 kV

### Max. Output Current

1 mA

### Max. Ch. Output Power

6 W

### Vset Resolution

20 mV

### Vmon Resolution

20 mV

### Iset Resolution

5 nA

### Imon Resolution

5 nA (high range) / 500pA (low range)

### Vmax

0 ÷ 6 kV

### Vmax resolution

2 V

### IMAX hardware

1 mA

### IMAX hardware resolution

1µA

### Ramp Up/Down

1 ÷ 500 Volt/s, 1 Volt/s step

## Trip

Max. time an "overcurrent" can last (seconds). A channel in "overcurrent" works as a current generator; output voltage varies in order to keep the output current lower than the programmed value. "Overcurrent" lasting more than set value (1 to 9999) causes the channel to "trip". Output voltage will drop to zero either at the Ramp-down rate or at the fastest available rate, depending on Power Down setting; in both cases the channel is put in the off state. If trip= INFINITE, "overcurrent" lasts indefinitely. TRIP range: 0 ÷ 999.9s; 1000 s = Infinite.  
Step = 0.1 s

## Voltage Ripple

10 - 1000Hz: <4mVpp typ; <8mVpp max  
1kHz - 20MHz: <2mVpp typ; <5mVpp max

## Vmon vs. Vout Accuracy

± 0.02% ± 1V

## Vset vs. Vout Accuracy

± 0.02% ± 1V

## Imon vs. Iout Accuracy

± 0.2% ± 1µA

## Iset vs. Iout Accuracy

± 0.2% ± 1µA

## Humidity range

0 ÷ 80% non condensing

## Operating temperature

0 ÷ 45°C

## Storage temperature

-10 ÷ 70°C

## Vout / Temperature coefficient

±50 ppm / °C

## Long Term stability (1 week after 1h warmup)

±0.02% / full scale

## Local Control

LCD touchscreen

## Remote Control

USB & Ethernet

## Ordering Options

Code	Description	
WDT8034XAAA	DT8034N - 8CH Desktop Programmable HV Power Supply(-6KV 1mA) -SHV conn. - Common Gnd	RoHS
WDT8034XMAAA	DT8034M - 8CH Desktop Programmable HV Power Supply(4ch +6KV 1mA, 4ch -6KV 1mA)-SHV conn. C. Gnd	RoHS
WDT8034XPAAA	DT8034P - 8CH Desktop Programmable HV Power Supply(+6KV 1mA) -SHV conn. - Common Gnd	RoHS

## Accessories


### HV CABLES



High Voltage Cable Assemblies

---

### A148x

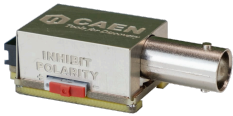

Inhibit - Kill Signal BNC Adapter for HV Power Supply Modules

---

### A1484

Inhibit - Kill Signal BNC Adapter for HV Power Supply Modules

---

### A1483

Inhibit - Kill Signal BNC Adapter for HV Power Supply Modules

---

## Related Software

### GECO2020



General Control Software for CAEN HV Power Supplies

---

## Related Software Libraries

### CAEN HV Wrapper Library



Library for CAEN Power Supply Control

## Related Products

### DT8033



8 Channel 4 kV/ 3 mA (6W) Desktop Power Supply Module (USB/Ethernet/Touchscreen)

### DT8031



8 Channel 100 V/ 10 mA Desktop Power Supply Module (USB/Ethernet/Touchscreen)

### DT8032



8 Channel 500 V/ 10 mA Desktop Power Supply Module (USB/Ethernet/Touchscreen)

### LabVIEW Driver (PSM - Power Supply Modules)



LabVIEW Instrument Driver for Power Supply Modules

**This document, or parts thereof, may not be reproduced in any form or by any means without written permission from Caen S.p.A. Although every effort has been made to ensure the accuracy of information presented in this catalog, Caen S.p.A reserves the right to modify its products specifications without giving any notice; for up to date information please visit [www.caen.it](http://www.caen.it) © Caen S.p.A - 2024**

**CAEN S.p.A.**

Via Vetraia 11  
55049 - Viareggio  
Italy

**Phone +39.0584.388.398**

**Fax +39.0584.388.959**

**info@caen.it**

**www.caen.it**

**CAEN GmbH**

Brunnenweg 9  
64331 Weiterstadt, Germany

**Phone +49 (0)212.254.4077**

**Mobile +49 (0)151.16.548.484**

**info@caen-de.com**

**www.caen-de.com**

**CAEN Technologies, Inc.**

1 Edgewater Street - Suite 101  
Staten Island, NY 10305  
USA

**Phone +1.718.981.0401**

**Fax +1.718.556.9185**

**info@caentechnologies.com**

**www.caentechnologies.com**

**CAENspa India Private Limited**

B205, BLDG42, B Wing,  
Azad Nagar Sangam CHS,  
Mhada Layout, Azad Nagar, Andheri West  
Mumbai, Maharashtra 400053, India

**info@caen-india.in**

**www.caen-india.in**

